

# Purple Vie® Thawer


## INNOVATION OF DRY THAWING SYSTEM

- Dry Thawing (Water-free)
- No-contact with direct heat
- AT-Closed vial only
- Reproducibility of result
- User friendly
- Traceability of data
- Portable (140 x 190 x 217 (mm))



## PERFORMANCE DATA

### ■ Water bath vs Purple vie

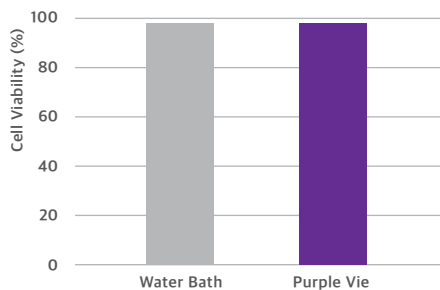


UCB-MS, CryoStor 10

### ■ Thawing Time (-80°C)

1.0mL	Approx 3min
1.5mL	Approx 3min 30sec
2.0mL	Approx 4min

AT-Closed vial 2.0mL, CryoStor 10

## HOW IT WORKS

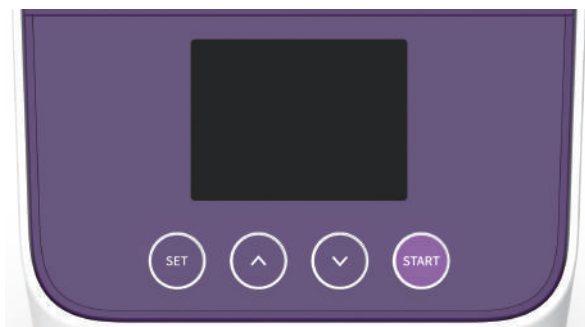

	SET	Setting of Thawing Time
	UP	10secs Up
	DOWN	10secs Down
	START	Thawing Start

## ORDERING INFORMATION

Product	Model. no.	Vial
Purple Vie® PV2 Thawer	PV2	2.0mL AT-Closed Vial
Purple Vie® PV6 Thawer	PV6	6.0mL AT-Closed Vial
Purple Vie® PV10 Thawer	PV10	10.0mL AT-Closed Vial

# *Purple Vie*

## **Manufacturer**

AMOLIFESCENCE Co., Ltd.

## **Technology Development**

AMOLIFESCENCE Co., Ltd., KANGSTEM BIOTECH Co., Ltd.

Website : <http://www.kangstem.com>

E-mail : [sales@kangstem.com](mailto:sales@kangstem.com)